



## ALCIAN BLUE STAIN pH 1.0 PROCEDURE

Item# STABL1

(Revised 04/28/11)

PRINCIPLE: This stain demonstrates strongly sulfated mucosubstances.

SPECIMEN: Any well fixed paraffin embedded tissue cut at 4 to 5 microns.

QUALITY CONTROL: American MasterTech Recommended Control Slides:  
Colon, CSC1025P; Small Intestine, CSS0325P

SOLUTION:

### **3% ACETIC ACID:**


Glacial Acetic Acid ..... 3 ml.  
Distilled Water ..... 97 ml.  
(Mix components *thoroughly* before use.)

PROCEDURE:

1. Deparaffinize slide using Xylene or a Xylene Substitute and hydrate through alcohols.
2. Rinse slide in running Tap water.
3. Rinse slide in Distilled water.
4. Place slide in **3% Acetic Acid** for 3 minutes.
5. Without rinsing, place slide in **Alcian Blue Stain** for 30 minutes.
6. Rinse slide quickly in running Tap water.
7. Place slide in **Nuclear Fast Red Stain** for 5 minutes.
8. Rinse slide in running Tap water for 1 minute.
9. Dehydrate slide through 3 changes of Absolute Alcohol.
10. Clear slide through 3 changes of Xylene or Xylene Substitute.
11. Coverslip using a permanent mounting media.

RESULTS:

Strongly sulfated mucosubstances: **BLUE**  
Nuclei: **PINK TO RED**  
Cytoplasm & Background: **PALE PINK**

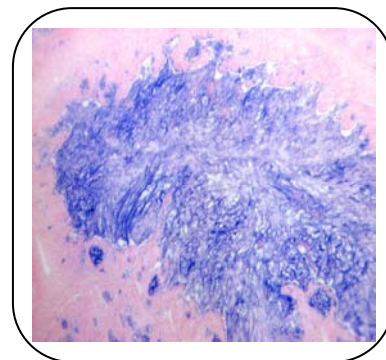

Human Colon Tumor

**REFERENCE**: Sheehan DC Hrapchak BB: Theory and Practice of Histotechnology; 1980; 172 - 173.

**Copyright ©1983-2012 American MasterTech Scientific Incorporated. All rights reserved.**

No part of these pages may be used for any purpose other than personal use. Therefore, reproduction, modification, storage in a retrieval system or retransmission, in any form or by any means, for reasons other than personal use, is strictly prohibited without prior written permission.